



Instructions to install the Image-Pro Plus driver for ProgRes® microscope cameras


1. Image Pro Plus 5.1 or higher must be previously installed.
2. Download the latest version of the Jenoptik Image-Pro driver for ProgRes® microscope cameras at our web site:
http://www.progres-camera.com/microscope-camera/downloads_software.html
3. Start the installation of Jenoptik Image-Pro driver for ProgRes® microscope cameras via double click on the icon and follow the instructions

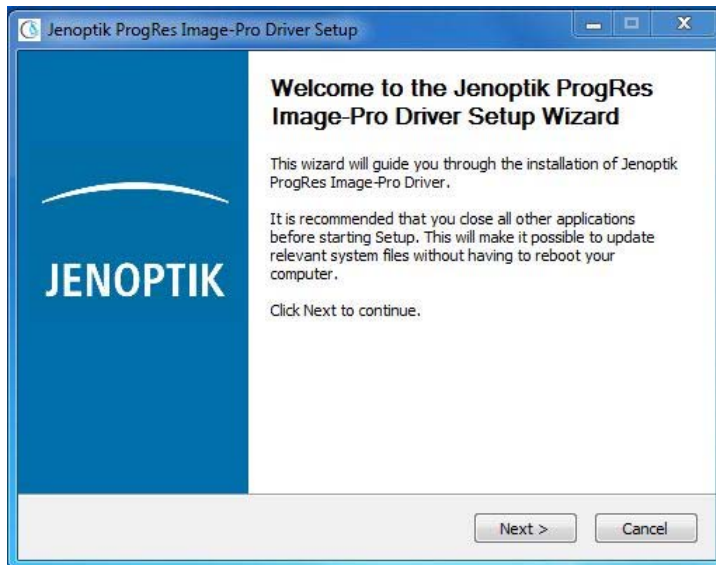


4. Please confirm the license agreement.

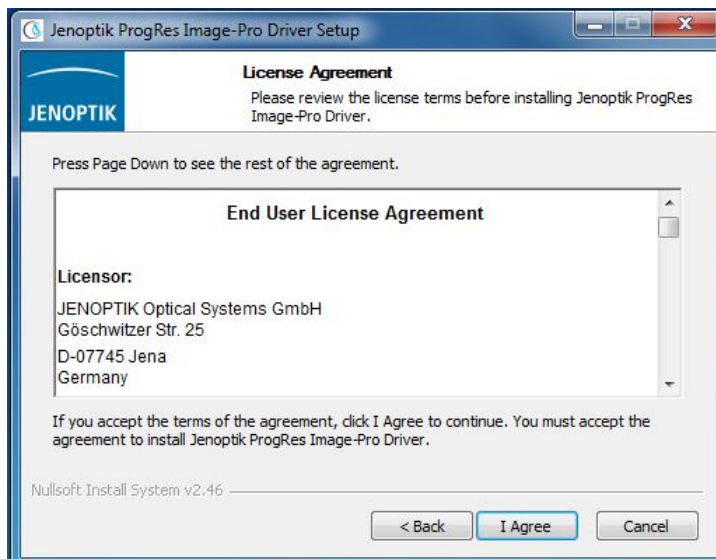

5. Please select the components to install.

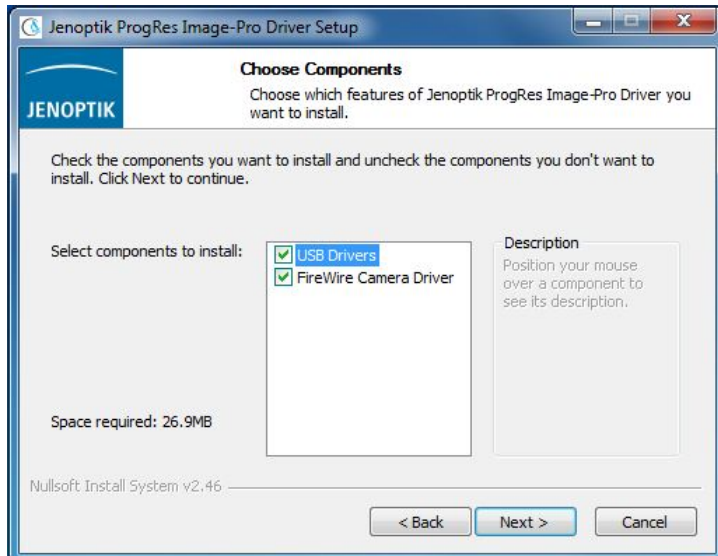


6. Please choose the Start Menu Folder.

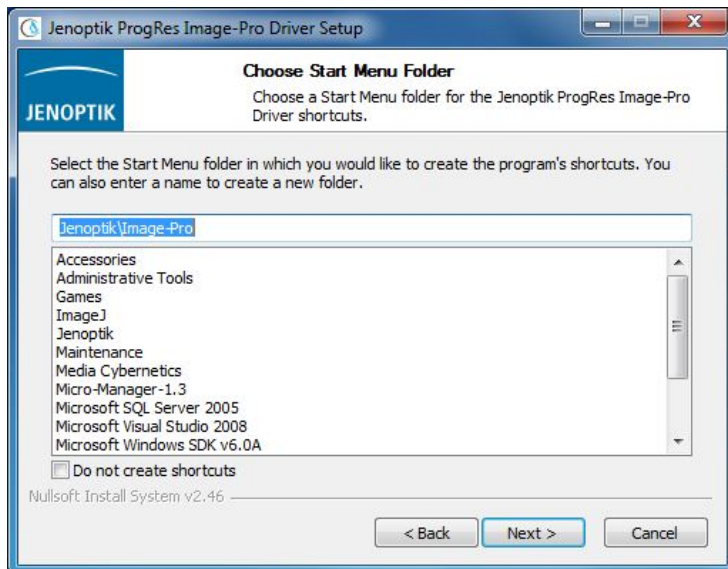


7. Please select the destination folder to install the Jenoptik Image-Pro driver for ProgRes® microscope cameras


Note: It is recommended to install the Jenoptik Image-Pro driver for ProgRes® microscope cameras in the Main folder of the Image-Pro Plus software.

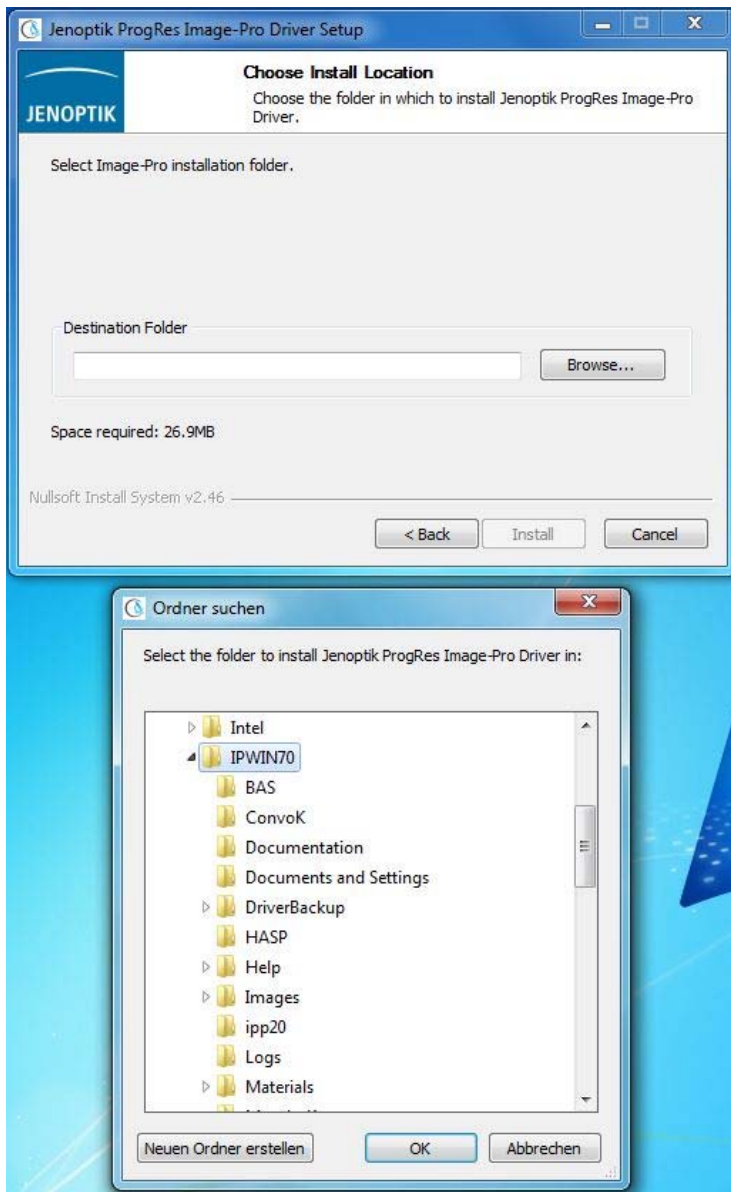

8. After the selection of the destination folder, please follow the instructions for installation.

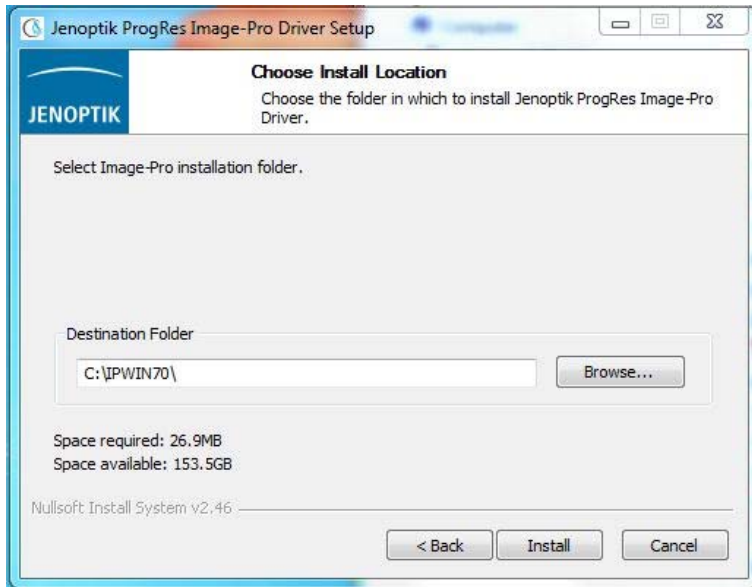


9. installation has been completed

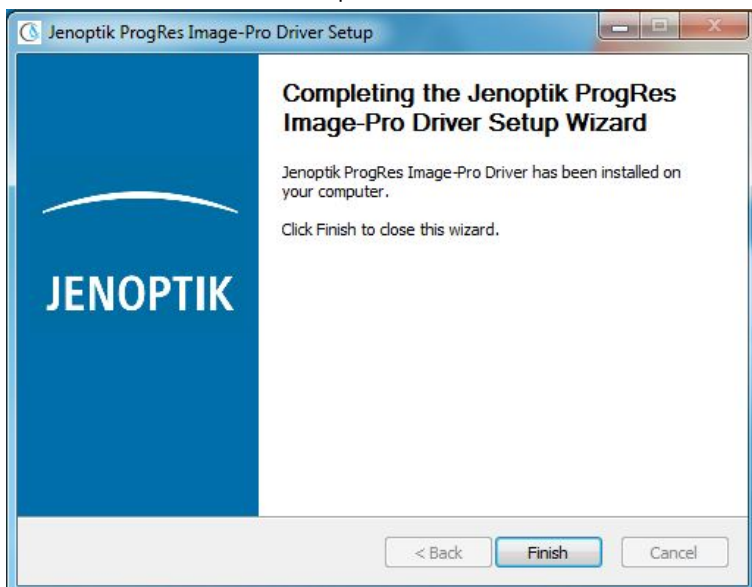

After connecting a ProgRes® microscope camera of Jenoptik, you are ready to run this camera with the Image-Pro Plus software.

10. Start the Image-Pro Plus software and the ProgRes® camera connection.

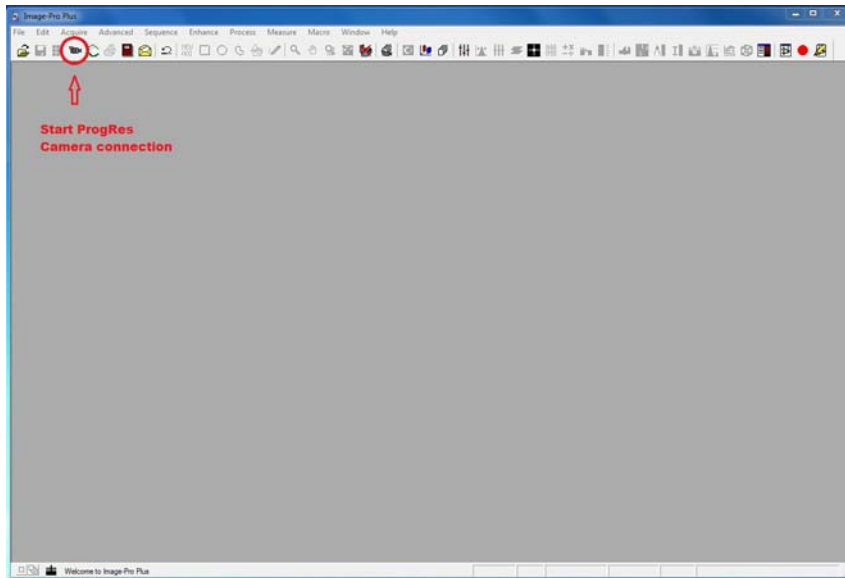

11. Choose the installed Jenoptik Image-Pro Plus driver.

