



Instructions to install the Jenoptik DirectShow driver for ProgRes® cameras

1. You can download the latest version of the Jenoptik DirectShow driver at our download area:
http://www.progres-camera.com/microscope-camera/downloads_software.html
2. **Attention: Please unplug all ProgRes® camera before you start with the installation of the DirectShow driver.**
3. Start the installation of Jenoptik DirectShow driver via double click on the icon and follow the instructions with "Next".




4. Please confirm the End User License Agreement with „I Agree“.


Note: The installation will be discontinued, if you disagree the End User License Agreement!



5. Please choose the components to install and follow the installation with „Next“.

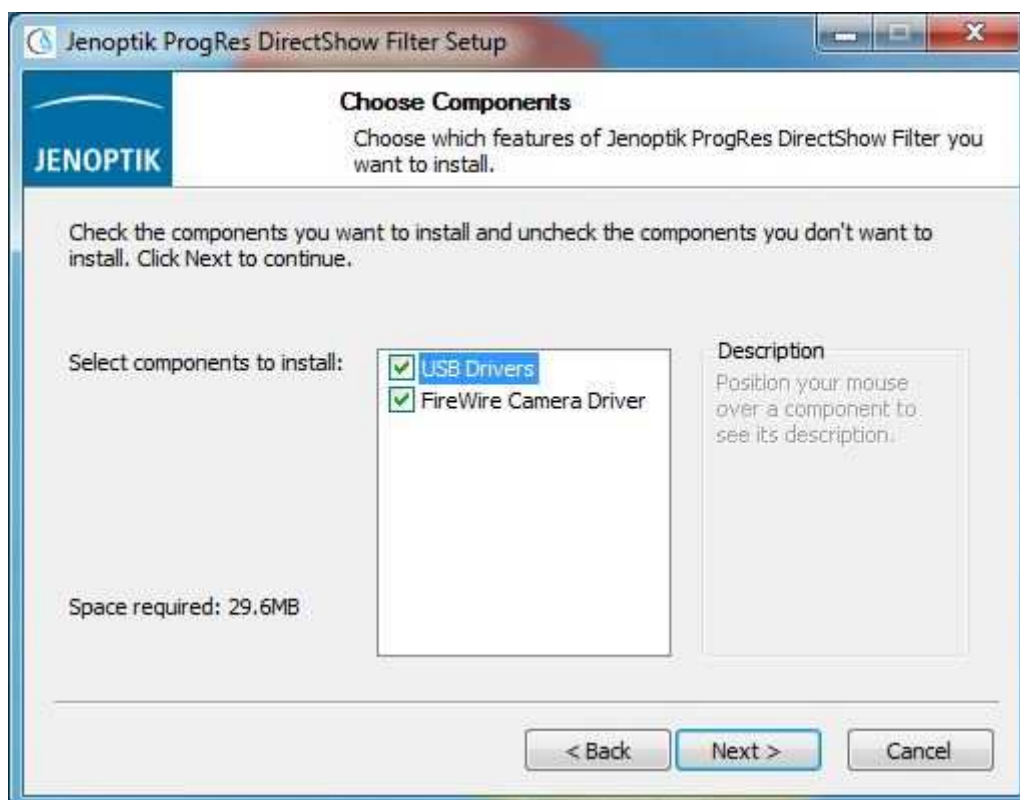

6. Please choose the install folder and follow the installation with „Next“.

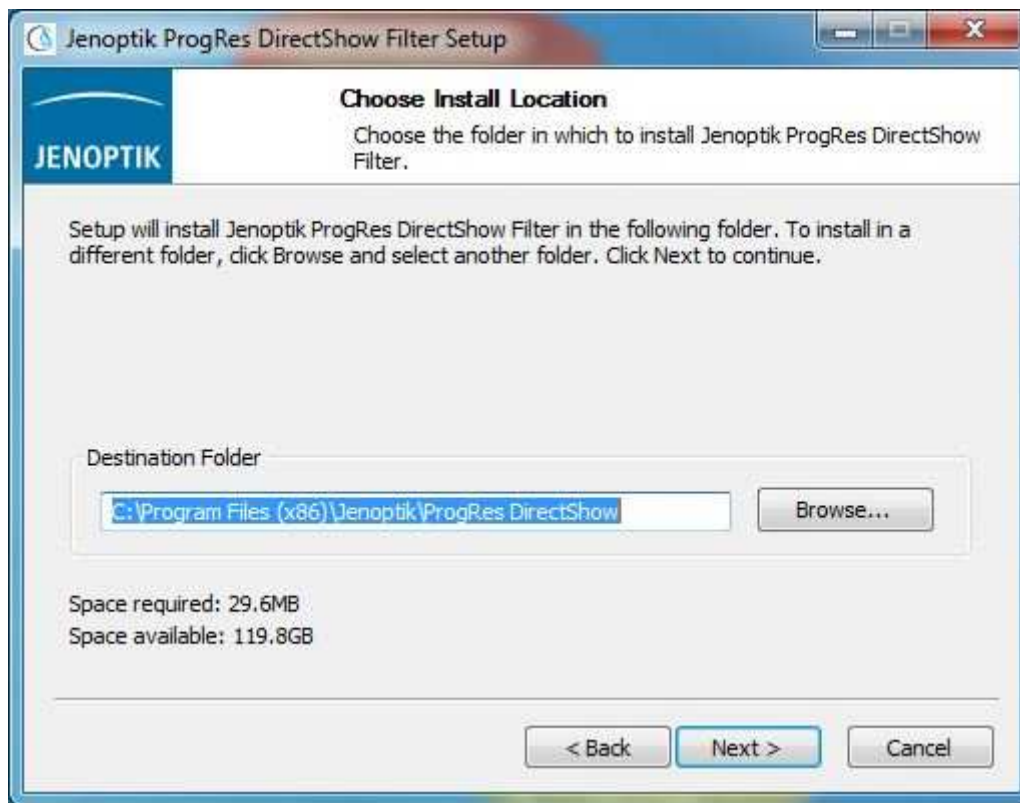


7. Please choose the Start Menu folder and follow the installation with „Next“.



8. If you want to use a ProgRes® camera with a Microsoft **64 bit** operating system, you have to install a special Firewire Bus driver previously.

Detailed information regarding the Firewire bus driver installation you can find under the following link:

<http://www.jenoptik.com/en-FAQ-DI-software-CapturePro-64-bit-busdriver>

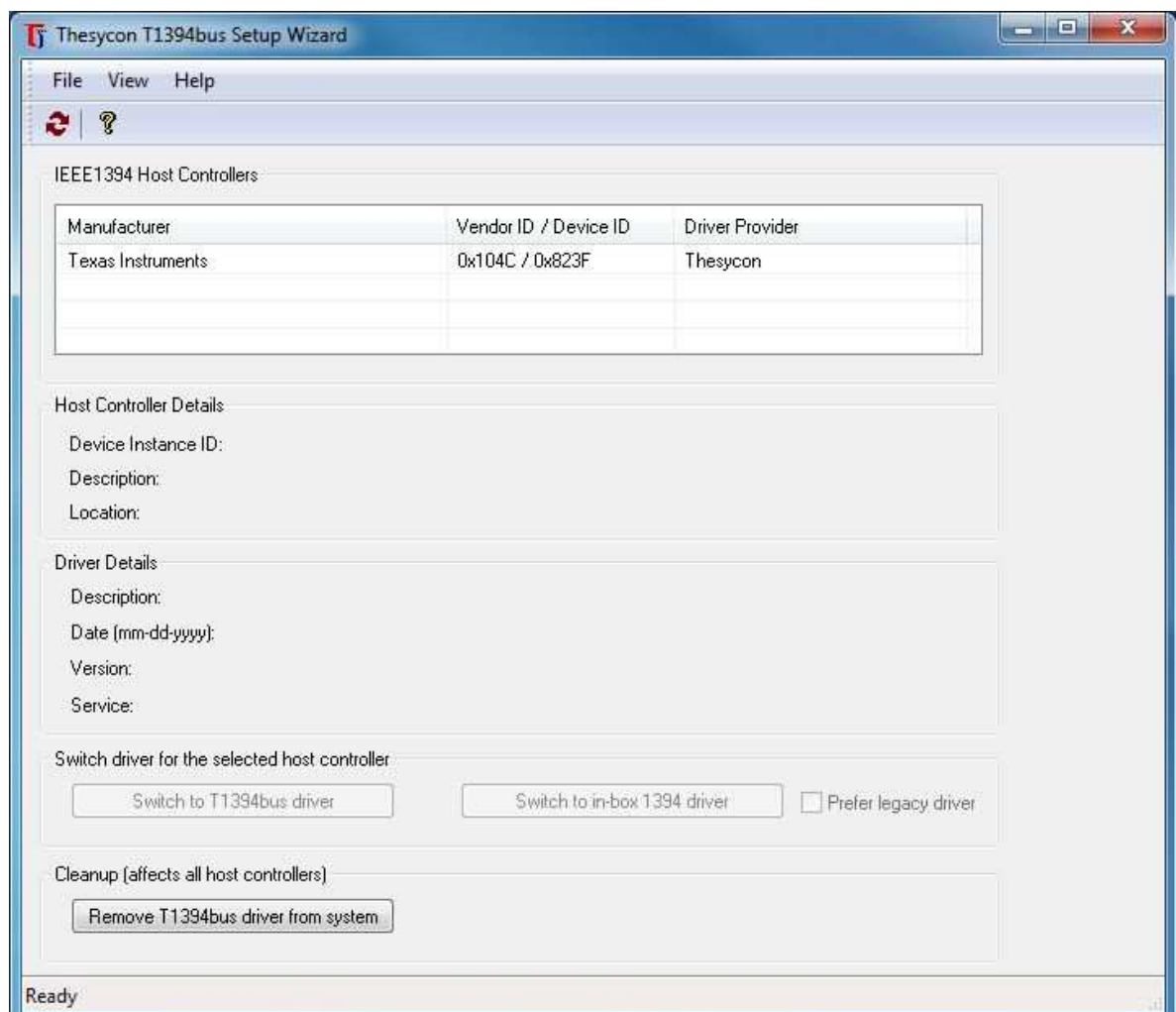

9. Please close the DirectShow installation after successfully installation of the camera device driver and bus drivers with "Finish".



10. Now you can plug in the ProgRes® camera to the computer and the Windows operating system will setup the driver automatically.
11. Note: You can find a list of supported cameras and resolution at the installation folder of the Jenoptik DirectShow driver.